



## Water Sampling Procedures

### WHIRL-PAK



1. Label the sample bag prior to filling. Permanent markers work great.
1. Fill the sample bag beyond the 4oz. fill line.
2. Pinch the top of the bag to close the gap.
3. Pinch the bag between the 100ml and 4oz. fill lines and pull gently to spill out the excess water.
4. Roll the wire tab towards the front of the bag until the bag is taut.
5. Wrap the wire tabs towards the back of the bag and twist to secure.
6. This procedure will leave the bag without any air in the bag.
7. Packaging and shipping
  - a. Pack all sample bags into a zip style plastic bag.
  - b. This bag should be packed gently inside the box with packaging material to prevent the samples from moving during shipment.

